

Installation

1. When unpacking, make sure the eyewash is intact.
Complete **KN 1724P**:
Eye wash, emergency sign and installation directions.
2. **The eye wash must have a flow of 28 litres per minute to function properly. The pipe dimensions must be adjusted to this flow!**

Water supply connection: ½" NPT female

Do not install the eye wash if correct flow can not be guaranteed!

The flow and pressure regulator of the eye wash work automatically at 2,4-6,5 bar. If the pressure exceeds 7 bar it will be necessary to install a reducing valve.
3. **Flush through the system thoroughly so that grindings and other particles are removed.**
4. Install the eye wash approximately 1100 mm from floor to spray heads.
5. Attach the enclosed emergency sign clearly visible.
6. Connect a waste pipe if necessary.
7. Check the function of the eyewash by flushing several times.
8. Hand over the instruction sheet to the safety supervisor or other responsible personnel.

Maintenance and operating instructions

1. Inform and train the personnel in question about the function of the eye wash and where it is located.
2. The eye wash is always in working order.

The eye wash activates by the push plate. The valve is not self closing, so the wash is turned off by pulling the plate back again.
3. The eye wash must be easily accessible for immediate use by the victim. Make sure that the area in front is not blocked.
4. **Practice the use of the device continuously.** By doing so, you will avoid stagnant water in the pipes. **Old water accumulates bacteria.**
5. **Use safety goggles.**